



PROCEDURE FOR COLLECTING A POOL SAMPLE

Note: Bottles are sterilized, do not touch the inside of the lid or lip of the bottle. Never open the bottle until collecting the sample. The powder or liquid that is inside the bottle is not to be rinsed out before sampling. This is a de-chlorinizing agent that is added for testing.

1. As far from the edge of the pool as possible, submerge bottle.
2. Once under water remove lid and fill bottle completely to the fill mark.
3. Replace lid while bottle is still under the water.
4. Take chlorine tests for free and total chlorine-record readings on data sheet. (This is optional and for your information only.)
5. Fill data tags out completely.

SPA SAMPLING

Use the same procedure as above when collecting your spa sample. Be sure to identify what location (ex. Pool or Spa) when filling out the data tags. A pseudomonas sample should also be collected from the spa. Pseudomonas is another type of bacteria which is common in spas. Once again, use the same sampling procedure as above.

1532 DeWitt Street
Ellsworth, IA 50075
Phone: 515 836.4444
Fax: 515 836.4541
lindaman@netins.net